

Adjusting Rinse Time



Compatible with software version 1.19.7 and newer



Please note:

- When idling, the instrument will automatically rinse the tubes with ddH₂O after a defined amount of time (**Rinse Time**). This prevents wash buffer crystallization.
- The default **Rinse Time** is 2 hours, but you can customize this setting.
- The instrument will rinse only **once** if no further action is taken.
- If you prime the tubes after rinsing, the instrument will rinse the tubes again after idling for the defined time period.

IntelliPlex 1000

Adjusting Rinse Time_P-6_V01


- Select **Setting**
- Select **Rinse Time**

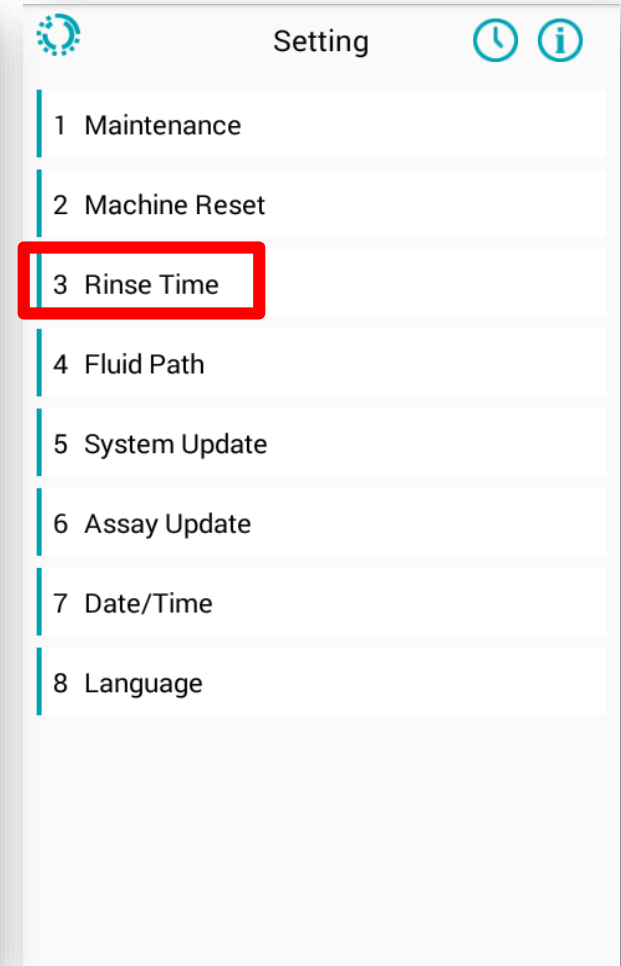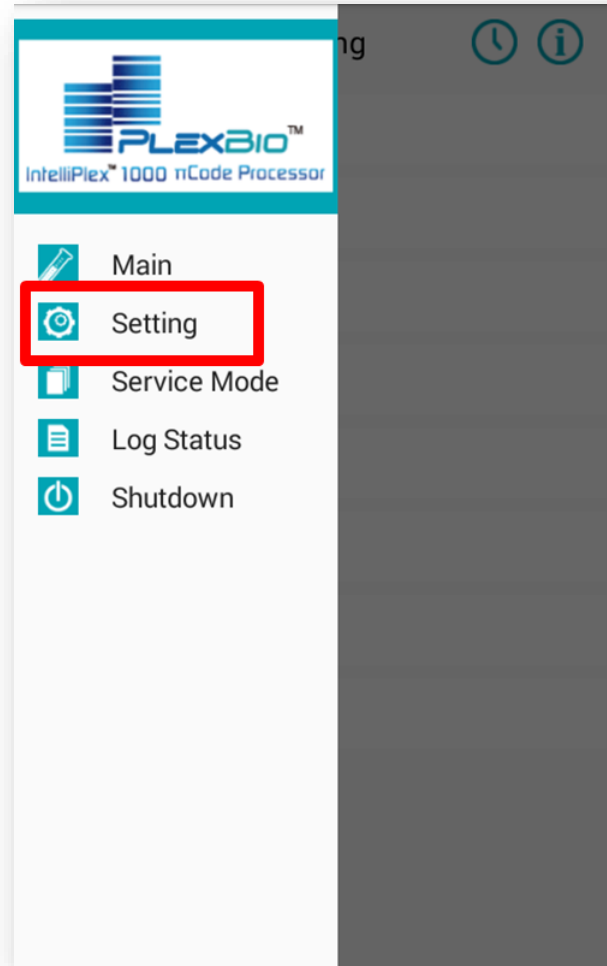


IntelliPlex 1000

Adjusting Rinse Time_P-6_V01

- Set the **Rinse Time** to the desired time interval (hh:mm:ss)

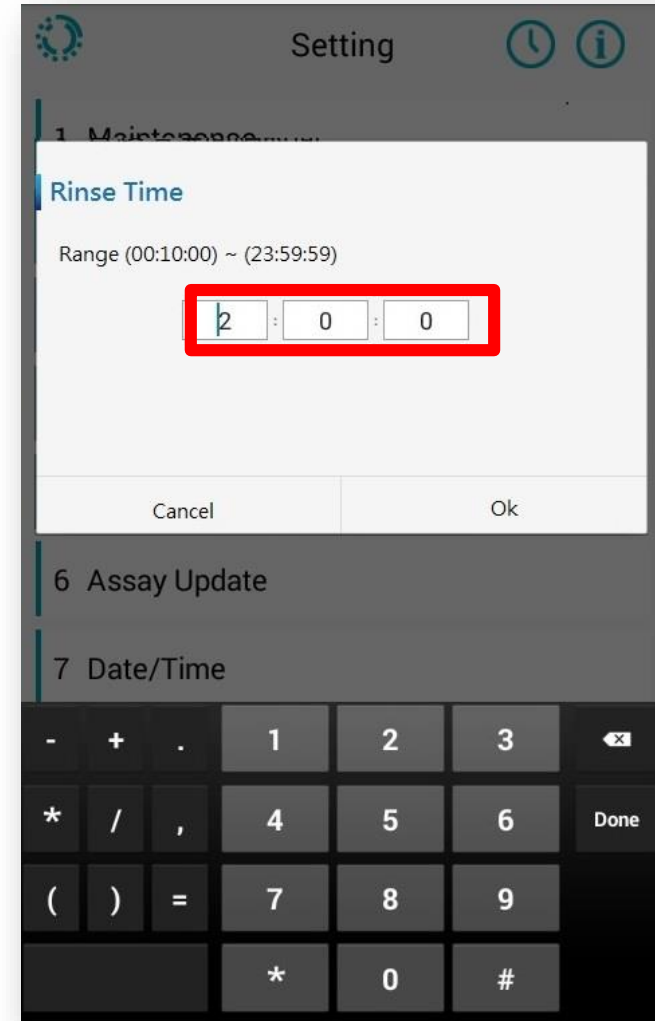

- Select **Ok**

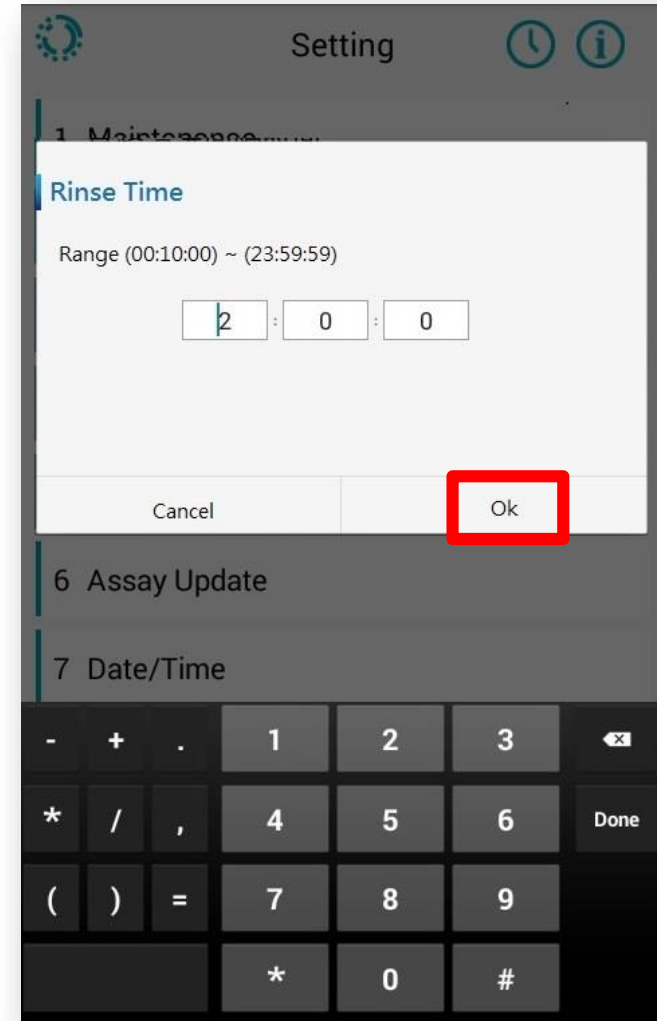

Thank You

PlexBio

6F-1, No. 351, Yangguang St.
Neihu District, Taipei City 11491, Taiwan
+866-2-2627-5878 | PlexBio.com
service@plexbio.com