

Exporting Raw Images


Software version 1.1.8.1429



Log in to DeXipher



Select Settings

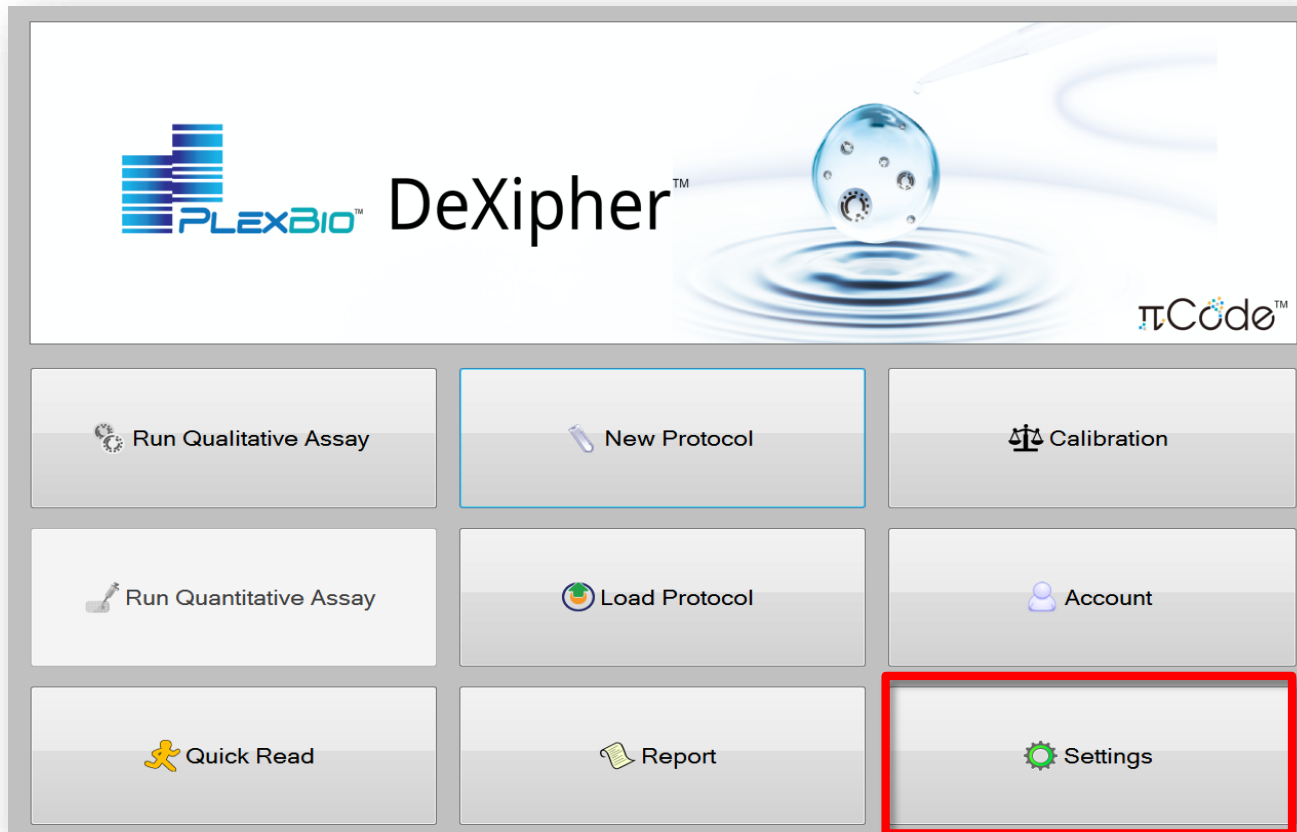


- Check the image file location


The screenshot displays the 'Report Setting' page in the PlexBio 100 software. On the left is a navigation sidebar with icons for Home, Quick Read, Run, Report, Calibration, Settings (highlighted), Account, and Logout. The main content area is divided into several sections:

- Report Setting:** Includes input fields for 'Min MicroDisc Count per πCode' (10), 'Max Replicate MFI CV' (100%), and 'Recovery Range' (± 30%). It also features dropdown menus for 'Standard Curve' (Best Fit) and 'Axis Transformation' (Log(x)-Linear(y)).
- Bar Chart Type:** Offers radio button options for 'MFI' (selected) and 'Concentration'.
- Report Columns:** A list of checkboxes for data columns: MFI - Bkgd, MFI SD, MFI CV, Dilution Factor, Obs. Conc., Unit, Exp. Conc., and (Obs/Exp)*100.
- MFI Section:** Contains checkboxes for 'MFI - Bkgd' (checked) and 'Show[MFI-Bkgd]<0 as 0'.
- Save Image:** A checkbox is checked and highlighted with a red box. Below it, the 'Image file location' is set to 'D:\plexbiolmage' with a browse button.
- Testing Lab Info:** Fields for 'Lab Name' and 'Lab Address'.
- View System Log:** A button with a circular arrow icon.
- Language Select:** A dropdown menu currently set to 'English' and a 'Save' button.

- Find the image file

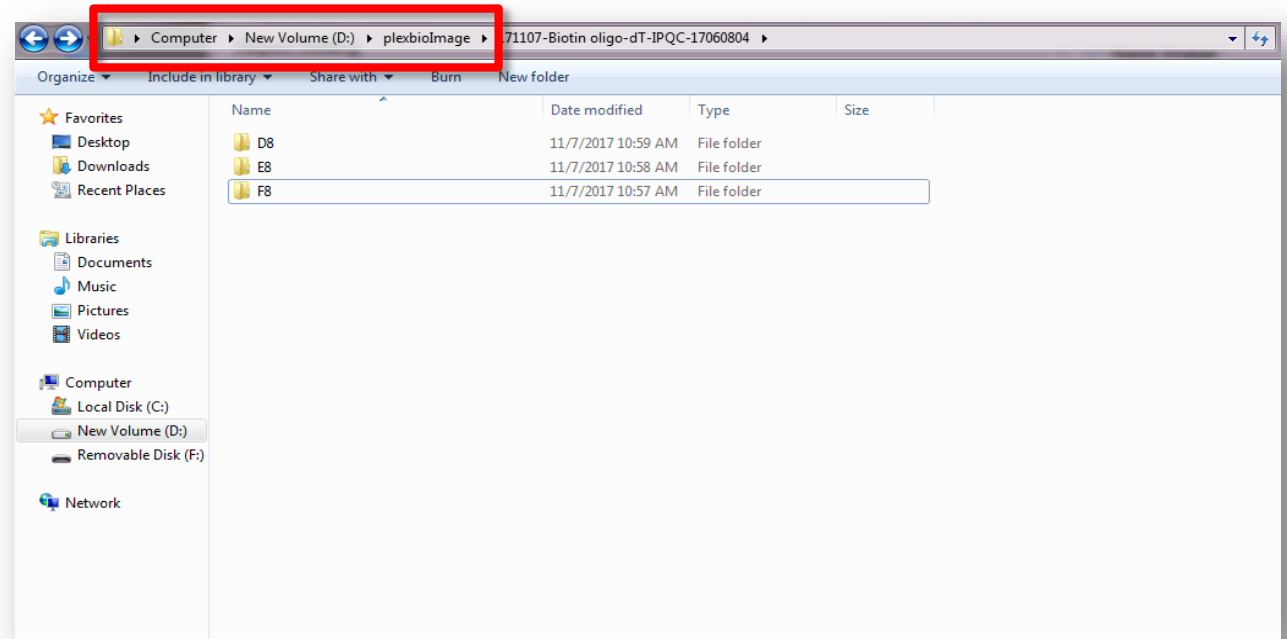

- Locate the image

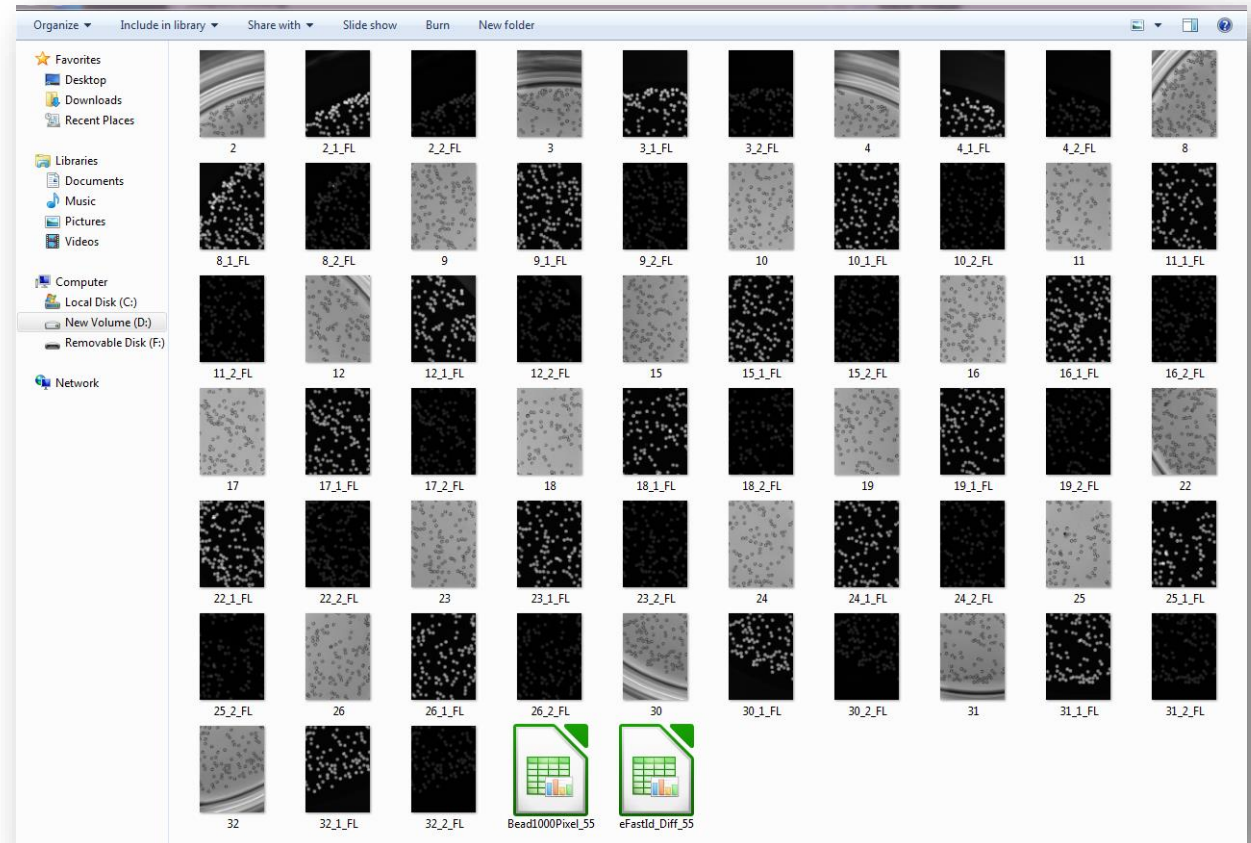

Thank You

PlexBio

6F-1, No. 351, Yangguang St.
Neihu District, Taipei City 11491, Taiwan
+866-2-2627-5878 | PlexBio.com
service@plexbio.com