

## Changing the DeXipher License

Software version 1.1.8.1429




## DeXipher Software License Features

Feature	Basic	RUO	MD
Account Management		✓	✓
Quick Read	✓	✓	
Qualitative Assay		✓	✓
Protocol Customization		✓	
Kit Report Generating		✓	✓
Fluorescence Calibration	✓	✓	✓
Electronic Signature (21 CFR Part 11)			✓


## Log in to DeXipher





- Select “i” for system information

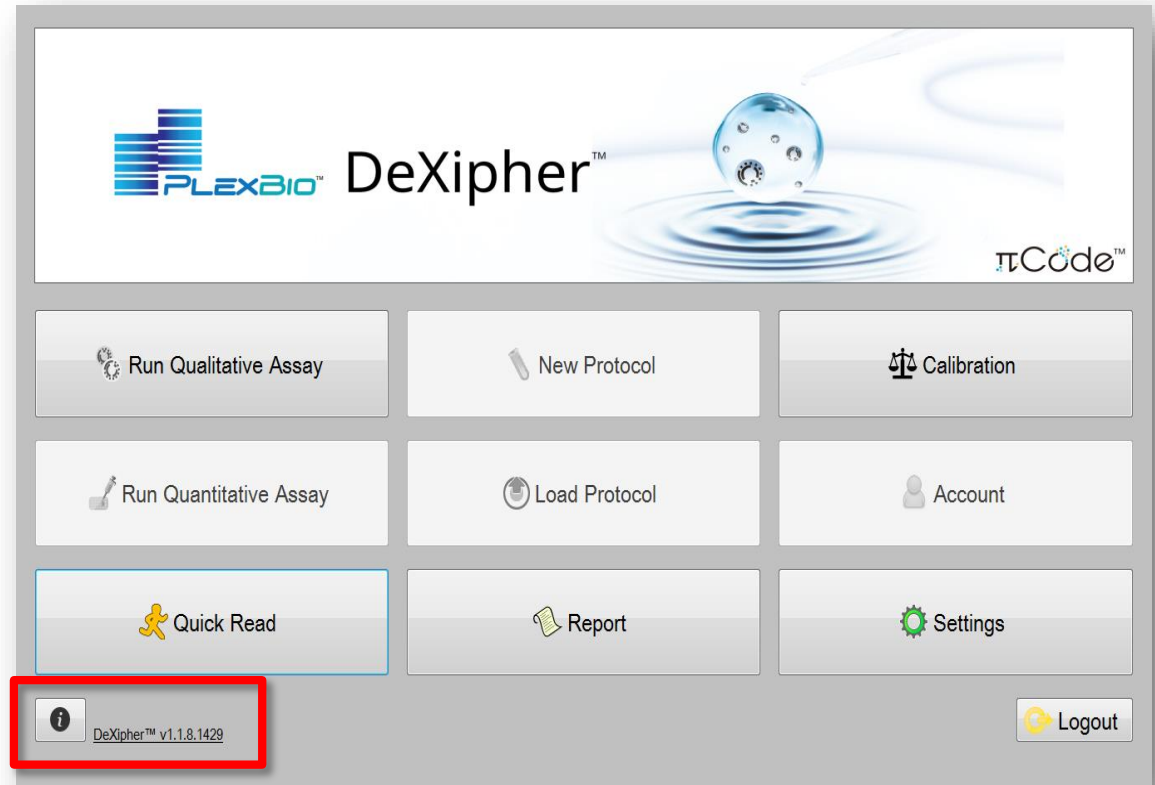


- Copy the Serial Number and send it to PlexBio.
- Receive the corresponding license key (purchased from PlexBio)

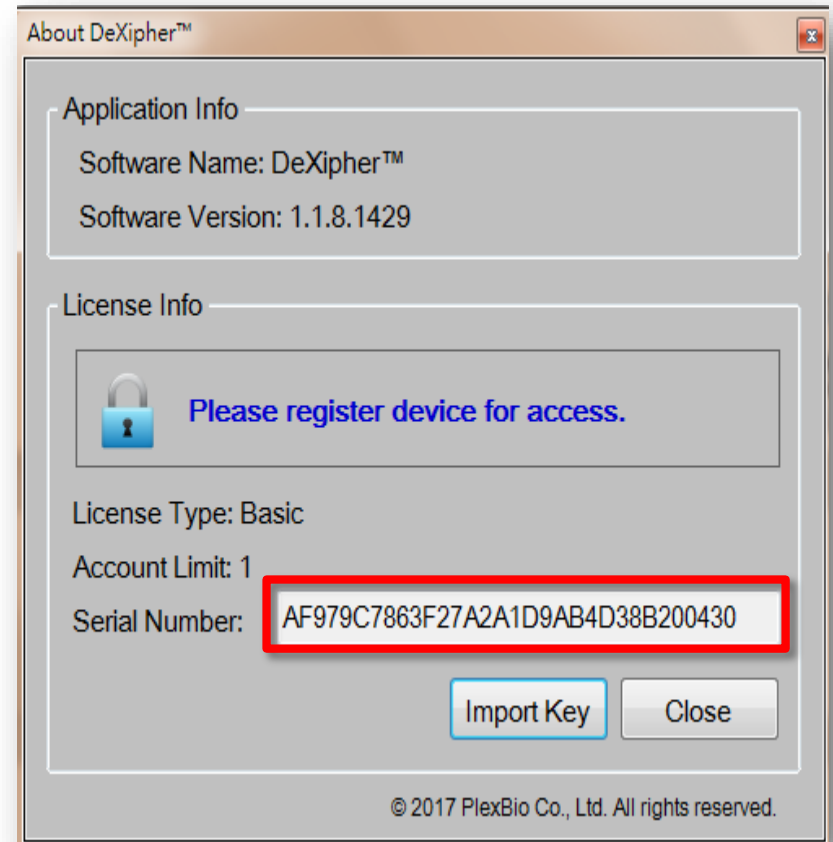


- Click **Import Key**
- Select the corresponding license key (acquired from PlexBio)

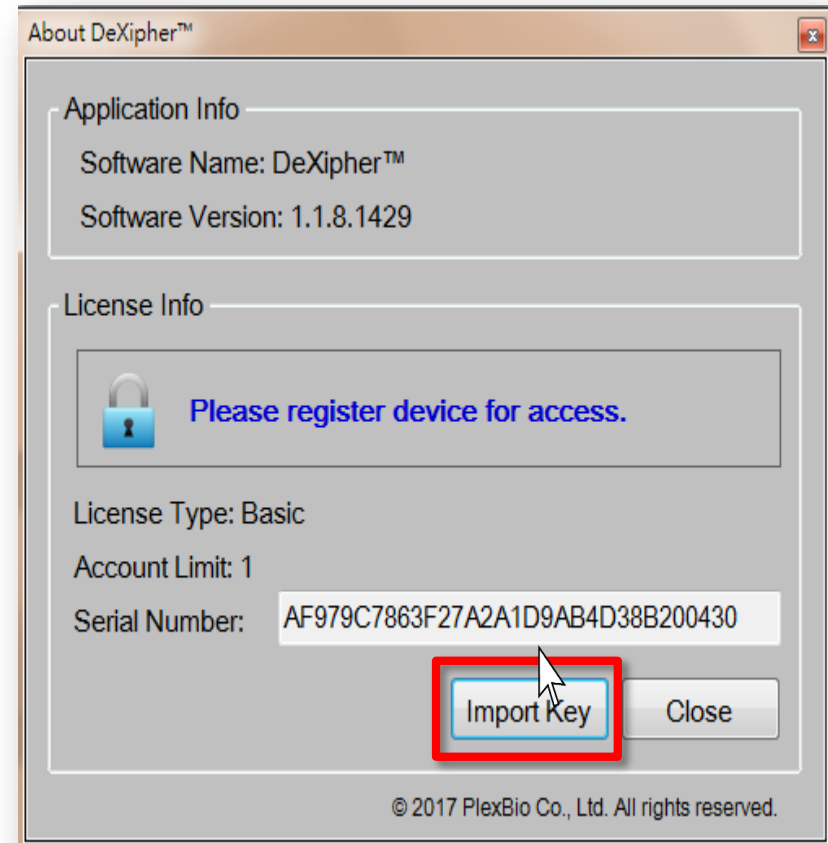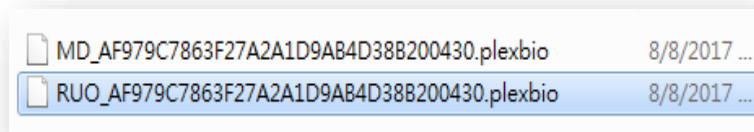

- Logout and login again

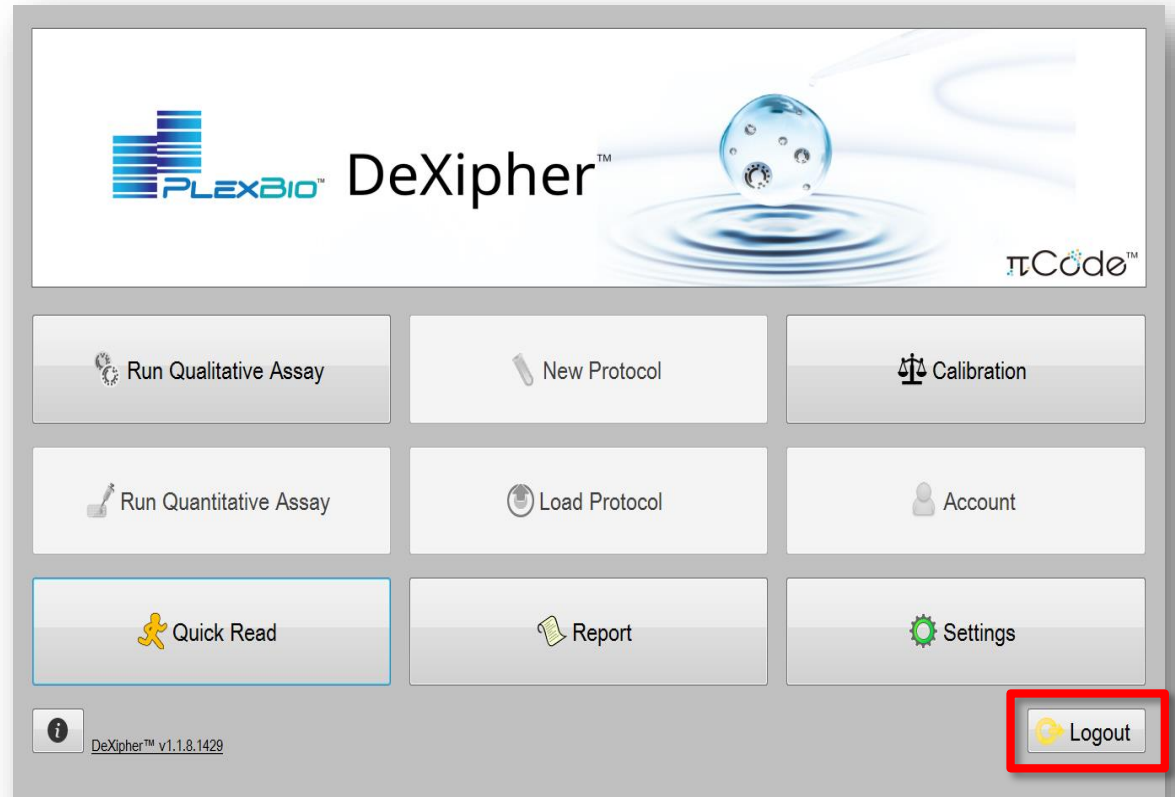

- Click **Information** and verify that the DeXipher license has been changed

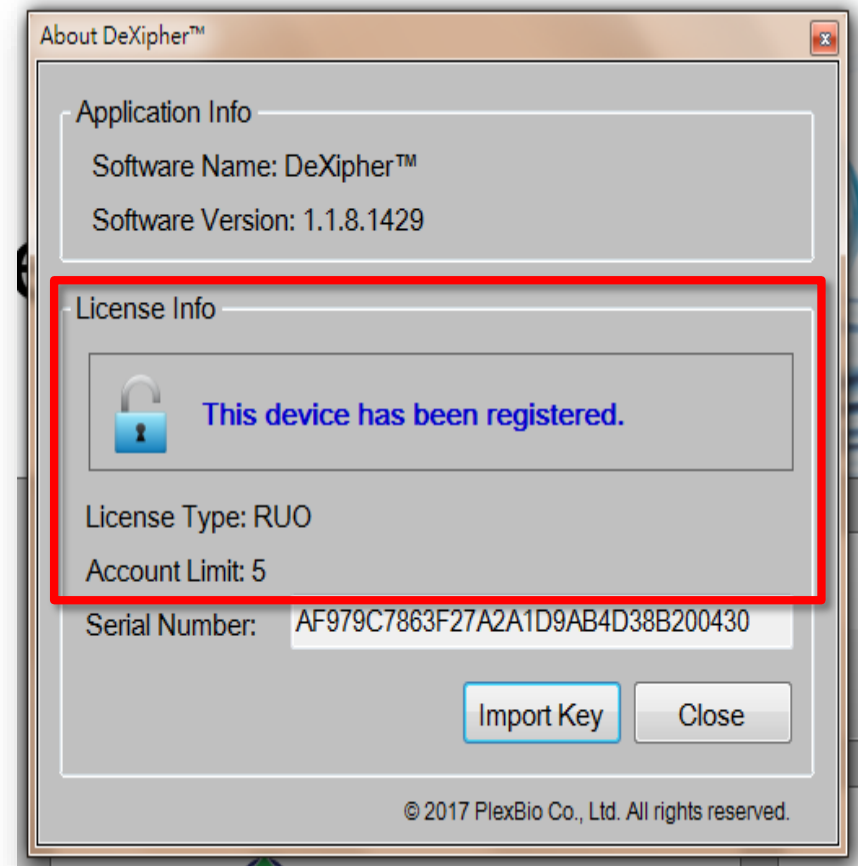



# Thank You

## **PlexBio**

6F-1, No. 351, Yangguang St.  
Neihu District, Taipei City 11491, Taiwan  
+866-2-2627-5878 | [PlexBio.com](http://PlexBio.com)  
[service@plexbio.com](mailto:service@plexbio.com)