

Importing an ENC File

Software version 1.1.9.1796


For RUO and MD license types only



Before you begin

- Ensure that the assay's Kit App has been installed
 - See **A-5_Installing Assay Kit Apps** for details

- Download ENC file
- Go to **[plexbio.com](https://www.plexbio.com)** and download the ENC file with the lot number that corresponds to your kit

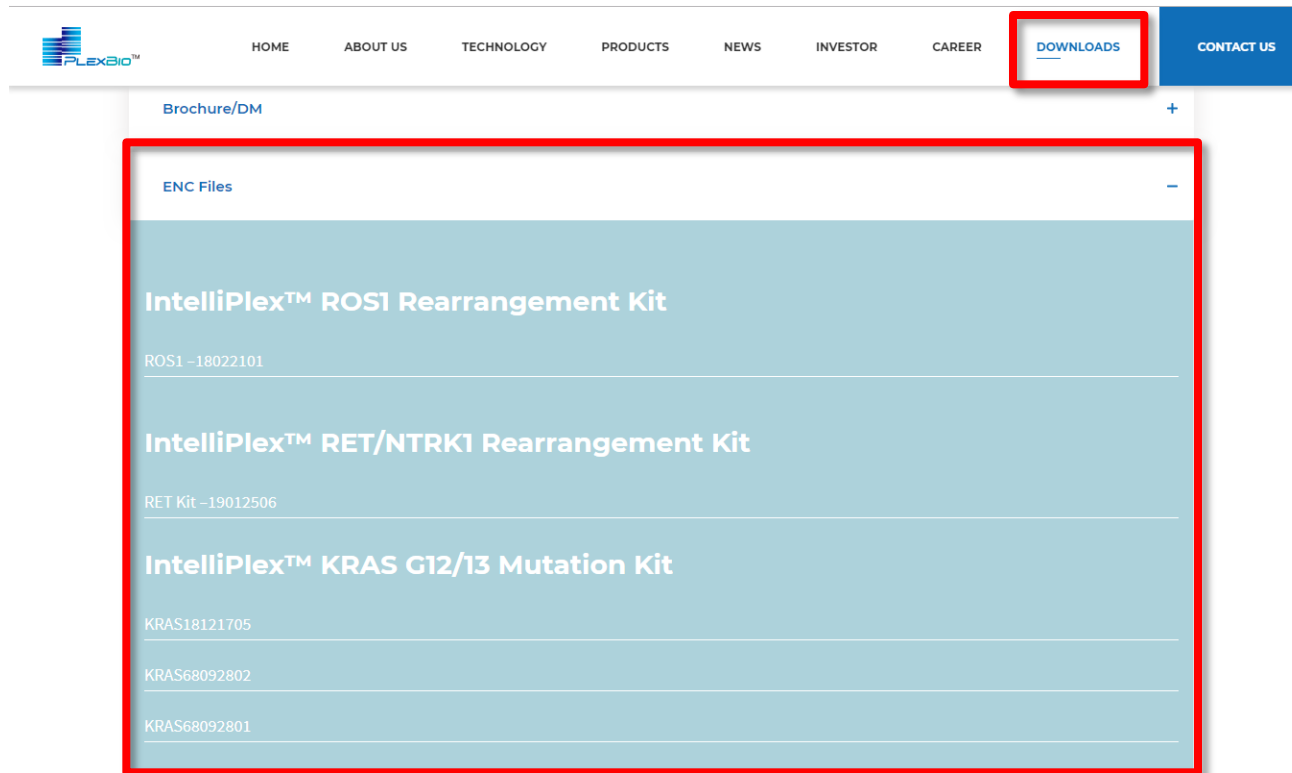

The screenshot shows the PlexBio website navigation bar with the 'DOWNLOADS' menu item highlighted in a red box. Below the navigation bar, a dropdown menu is open, showing a list of ENC files. The 'ENC Files' section is also highlighted in a red box. The list includes:

- IntelliPlex™ ROS1 Rearrangement Kit
ROS1 -18022101
- IntelliPlex™ RET/NTRK1 Rearrangement Kit
RET Kit -19012506
- IntelliPlex™ KRAS G12/13 Mutation Kit
KRAS18121705
KRAS68092802
KRAS68092801


PlexBio 100

Importing an ENC File_A-6_V02


- Log in to DeXipher



- Click **Run Qualitative Assay** or **Run Quantitative Assay**

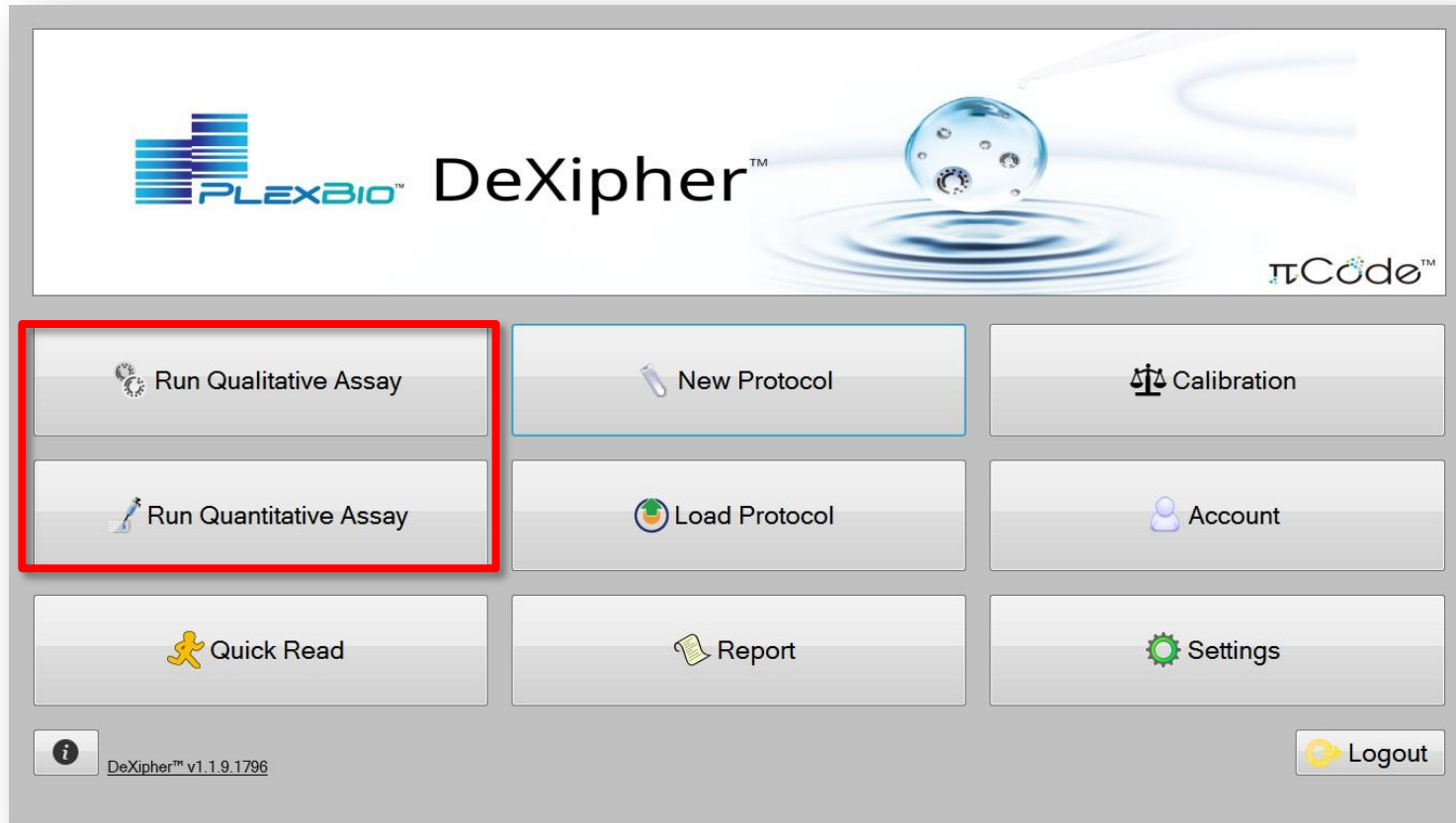

- Click **Import Kit** and Select the corresponding ENC file to import

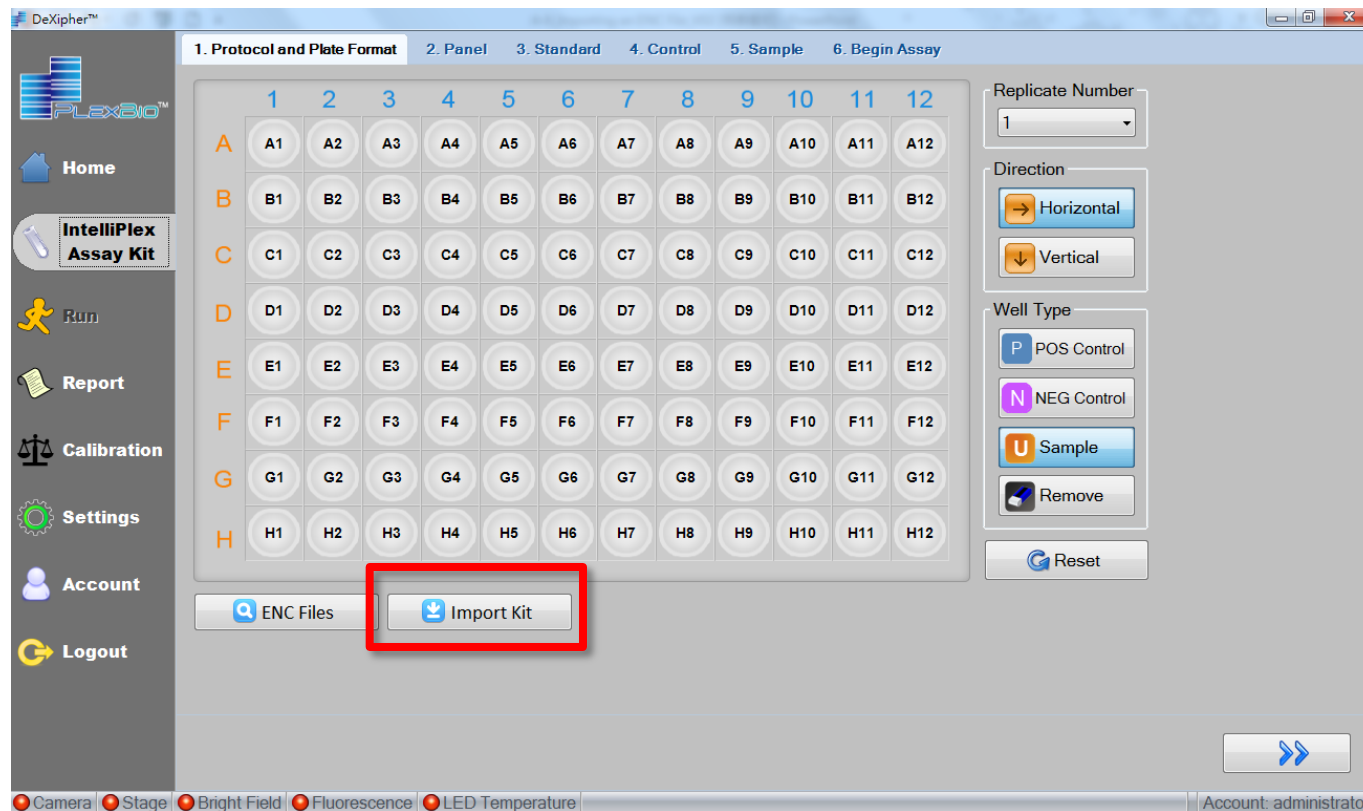

- Click **ENC Files** to check if the corresponding ENC file has been imported.

The screenshot shows the DeXipher software interface. The main window displays a 96-well plate layout (A1-H12) and various assay settings. A dialog box titled 'ENC Files' is open, showing a table of imported kits. The 'ENC Files' button in the main interface is highlighted with a red box.

Kit Name	Lot No.	Expiry Date	Import Date
HCV Kit	18010501	2018/08/25	2015/04/24
HCV 16111602	16111602	2017/05/16	2017/02/15
HCV 16111601	16111602	2017/05/16	2017/02/24
HCV Genotyping Kit	17020601	2017/07/06	2017/04/10
KRAS17071104	17071104	2017/10/31	2017/07/12
KRAS17092201	17092201	2018/07/22	2017/09/25
KRAS17111001	17111001	2018/09/03	2017/11/13
KRAS17111703	17111703	2018/09/03	2017/11/20
APP HPV Kit ENC File	17081802	2017/10/18	2017/11/23
KRAS17112301	17112301	2018/09/21	2017/11/27
HCV Kit - 17080301	17080301	2019/12/31	2018/01/10
HCV Kit - 18010501	18010501	2018/08/25	2018/01/10

Thank You

PlexBio

6F-1, No. 351, Yangguang St.
Neihu District, Taipei City 11491, Taiwan
+866-2-2627-5878 | PlexBio.com
service@plexbio.com